Instrumentation

Alimentation

Instrumentation

Power supply

Ref :
281 084

Français – p 1

English – p 5

Version : 6009

Alimentation Evolution[®] F15 +/-
Evolution[®] power supply F15 +/-

ALIMENTATION STABILISEE - 15, 0, + 15 VOLTS / 0,5 AMPERE

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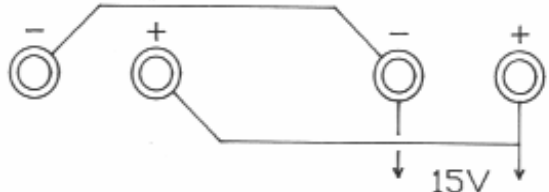
STABILISED SUPPLY - 15, 0, + 15 VOLTS / 0.5 AMP

1 Use

- Plug into the 230 volt mains supply.
- Put the general switch to "ON". The "mains on" indicator lights.
- Connect your circuit according to the use desired.

	2 x 15 V / 500 mA supplies
	- 15, + 15 V / 500 mA symmetrical supply
	30 V / 500 mA supply
	15 V / 1 A supply

PRECAUTIONS IN USE

- Never obstruct the ventilation apertures of the power supply when it is in use.
- Never use the power supply in the immediate proximity of liquids (water, acids, etc.).
- Avoid leaving the power supply running in full sun or near a significant heat source.

Keep the original packaging; it will be required to return the equipment to our factory for any repairs needed.

2 Technical features

Complies with the EN61010-1 norm.
 Category II, degree of pollution 2.

2.1 Input

- Mains supply 230 volts $\pm$ 10 %, 50-60 Hz.
- No earth terminal (double insulated).
- Consumption: 35 VA max.
- Protected by 315 mA slow fuse (not accessible to the user).

2.2 Outputs

- 2 independent 15 V, 500 mA floating outputs connectable in series or parallel according to use.
- Stability greater than 0.1 %.
- Residual ripple less than 5 mV.
- Protection by current limitation.

2.3 Mechanical features

- Stackable casing in shock resistant plastic (ABS high temperature, self extinguishing).
- Ambient operating temperature: 0 to 40 °C.
- Dimensions of casing: 230 x 230 x 125 mm.
- Weight: 1.8 kg.

3 Care and maintenance

In instances of non operation:

- Check that the mains supply is working (230 V).
- Check your test piece so as to detect any possible short circuit or flow of current greater than 500 mA (power supply tripped).

4 After-Sales Service

This material is under a two year warranty and should be returned to our stores in the event of any defects.

For any repairs, adjustments or spare parts, please contact:

JEULIN - TECHNICAL SUPPORT
Rue Jacques Monod
BP 1900
27 019 EVREUX CEDEX FRANCE
+33 (0)2 32 29 40 50

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Service gratuit * :
+ 33 (0)2 32 29 40 50

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Rubrique FAQ

Direct connection for technical support

A team of experts at your disposal from Monday to Friday (opening hours)

- You're looking for technical information ?
- You wish advice for use ?
- You need an urgent diagnosis ?

We take in charge your request immediately to provide you with the right answers regarding your activity field : Biology, Physics, Chemistry, Technology .

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