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Small equipment

**Ref :
703 172**

Français – p 1

English – p 3

Version : 6007

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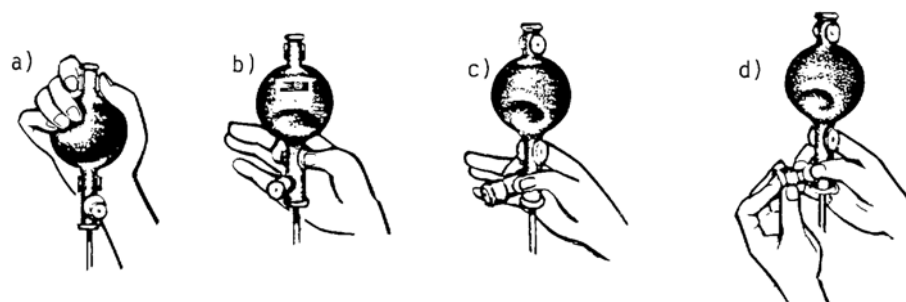
NOTES

The « ASSISTENT » pipette ball has 3 outlets, is produced without seams and therefore can be cleaned with acetone and other solvents.

1 Instruction for use

- a) The ball is held in the palm of the hand, thumb and forefinger on outlet R, and air is pressed out of the ball.
- b) Thumb and forefinger are pressed lightly on outlet A (indication valve) to draw in the required amount of liquid.
- c) For draining-off of liquid thumb and forefinger is held over outlet M.
- d) To achieve a complete and fast drainage, outlet M is pressed, the opening closed with the forefinger of the left hand, and in addition the olive is pressed between the thumb and middle finger

Note

In the case of liquid appearing in the ball, press out valve-ball "R" with a ruler-edge, then push out valve-ball "A" with a metal-rod of 4-5 mm diam. Wash and dry the pipette-ball, then replace valve-balls in their position. Avoid getting valve-balls inside of the pipette-ball.

2 After-sales service

The device is under a 2-year guarantee, it must be sent back to our workshops.

For any repairs, adjustments or spare parts please contact:

JEULIN – TECHNICAL SUPPORT
Rue Jacques Monod
BP 1900
27 019 EVREUX CEDEX FRANCE
+33 (0)2 32 29 40 50

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SA capital 3 233 762 € - Siren R.C.S. B 387 901 044 - Siret 387 901 04400017

Direct connection for technical support

A team of experts at your
disposal from Monday
to Friday (opening hours)

- You're looking for technical information ?
- You wish advice for use ?
- You need an urgent diagnosis ?

We take in charge your request immediatly to provide you with the right answers regarding your activity field : Biology, Physics, Chemistry, Technology .

Free service * :
+ 33 (0)2 32 29 40 50

** Call cost not included*