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Measurement and instrumentation

**Ref :
251 039**

Français – p 1

English – p 3

Version : 8010

Pressiomètre Initio®
Initio® pressure meter

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1 Technical specifications

- Range: 0 to 2000 hPa
- Accuracy: 2% +/- 4 hPa
- Resolution: 1 hPa
- Maximum pressure: 4000 hPa
- Dimensions: 100 x 100 x 40 mm
- Power supply: 1 6F22 type 9V battery

2 Precautions for use

Never introduce liquid or corrosive gas into the pressure sensor.

3 Replacing the battery

When the "Bat" indicator lamp comes on at the top left of the display, it is recommended that the battery be replaced.

To open the battery-housing, insert a small screwdriver in the notch at the edge of the flap and lever it up.

4 Offset adjustment

Adjusting the displayed value: Although it is set at the workshop, it is possible to modify the value displayed using the potentiometer located at the back of the battery compartment.

To do this, insert a small screwdriver in the RH hole and rotate till you obtain the desired value (for example, by keeping the sensor inlet in ambient air and by comparing the displayed value to the day's atmospheric pressure). This adjustment corresponds to an offset adjustment, and will offset all the measurements by more or less, by the same value.

5 After-sales service

The device is under a 2-year guarantee, it must be sent back to our workshops.

For any repairs, adjustments or spare parts please contact:

JEULIN – TECHNICAL SUPPORT
Rue Jacques Monod
BP 1900
27 019 EVREUX CEDEX FRANCE
0825 563 563

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We declare under our own responsibility, that the following products :

**INITIO ref : 302 174, 291 194, 291 195, 701 296, 251 039, 251 040, 251 041,
222 032, 222 035, 302 183, 302 184, 272 014, 272 015, 251 045, 701 332**

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H YASSINE

Responsable Technique

A stylized, handwritten signature in blue ink, consisting of several loops and a long horizontal stroke.

J-Y FRIBOULET

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A handwritten signature in blue ink, featuring a large, open loop at the beginning and a long horizontal stroke.

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SA capital 3 233 762 € - Siren R.C.S. B 387 901 044 - Siret 387 901 04400017

Direct connection for technical support

A team of experts at your
disposal from Monday
to Friday (opening hours)

- You're looking for technical information ?
- You wish advice for use ?
- You need an urgent diagnosis ?

We take in charge your request immediatly to provide you with the right answers regarding your activity field : Biology, Physics, Chemistry, Technology .

Free service *

+ 33 (0)2 32 29 40 50**

* Call cost not included

** Only for call from foreign countries